

## Setaflash™ Series 3

### The safest choice

Approved in over 1000 international product specifications and regulations

- ISO 3679; ISO 3680 (obs); ISO 9038
- ASTM D3278; D3828; D4206; D7236; EPA 1020 A & B
- IP 523; 524; 534; 602; BS 2000-523



1 minute test • 2ml sample • 0 to 300°C • Portable • Conforms



## Why use Setaflash Series 3?

- Conforms to international regulations
- 1 or 2 minutes test time
- 2ml sample
- 0°C to 300°C

## Key Features

- Compact, portable and rugged design
- Best published precision of any flash point method
- Suitable for unknown samples using ramp mode
- Automated flash detection
- Automatic barometric pressure correction
- USB port and result storage
- Low cost of purchase and maintenance



## Where can I find more Information?

Learn more about Setaflash testing:  
[www.stanhope-seta.co.uk/small-scale-flashpoint-testing.asp](http://www.stanhope-seta.co.uk/small-scale-flashpoint-testing.asp)



Watch the Series 3 video demonstration:  
[www.stanhope-seta.co.uk/5134/Setaflash-Series-3-Flash-Point-Tester](http://www.stanhope-seta.co.uk/5134/Setaflash-Series-3-Flash-Point-Tester)




## Operator Sequence

4 easy steps:



Select test temperature

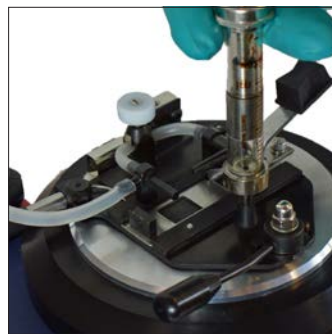

Inject 2ml sample



Press 




Dip the test flame, flash detection is automatic



## Trusted Precision & Accuracy

- The most precise flashpoint test
- Proven performance
- Approved in an extensive range of specifications and regulations

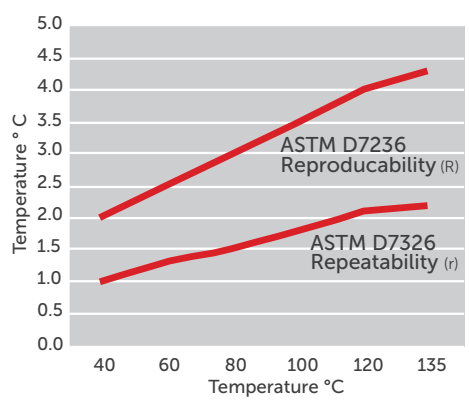

## USB Data Transfer

- 1GB data capacity, holding 100,000 results
- Results can be saved to a USB stick



## Calibration and Verification

- Software enables the user to calibrate temperature and barometric pressure
- Verification through Seta's Small Scale Certified Flash Point material (reference 99878-3)



## Small Sample - 2ml

- Small sample size of only 2ml
- Reduces costs and waste per test



## Service and Maintenance

- Rugged design
- Minimum maintenance
- Visit our technical support website: [www.stanhope-seta.co.uk/support](http://www.stanhope-seta.co.uk/support)



## Rapid Results in just 1 or 2 Minutes

- Fastest method of checking flashpoint
- Results determined in just one or two minutes
- Maximises operability



## Search for the Flash Point

- Test unknown samples using ramp mode
- Ramp mode is integral to 33000-2 model
- User adjustable ramp rates and dip interval

Technical Specifications:		
Seta Part Number:	30000-2	33000-2
Cup:	Closed	Closed
Material:	Aluminium	Aluminium
Temperature Range:	Ambient to 300°C	Ambient to 300°C
Temperature Ramp Mode:	N/A	Automatic or custom settings (ramp rate: up to 6°C/min and beep rate: up to 5°C/B intervals). Extended test range, 100°C from start temperature
Temperature Selection:	°C or °F	°C or °F
User Interface:	Colour LCD display - touch screen sensitive	Colour LCD display - touch screen sensitive
Navigation:	Control knob (rotational and push-down)	Control knob (rotational and push-down)
Data Storage and Download:	1GB internal memory and integrated USB port	1GB internal memory and integrated USB port
Fuel Supply:	Integrated gas tank (removable)	Integrated gas tank (removable)
Sample Size:	2ml or 4ml	2ml or 4ml
Pressure Correction:	Automatic barometric correction	Automatic barometric correction
Voltage Supply:	110 to 250V a.c.-50/60 Hz (universal)	110 to 250V a.c.-50/60 Hz (universal)
Power:	200W (Max.)	200W (Max.)
Size (LxWxH) / Weight:	28 x 14 x 21cm / 3Kg	28 x 14 x 21cm / 3Kg

## Products & Specifications

Diesel | Aviation Fuel | Marine Fuel | Biodiesel | FAME | Chemicals | Pharmaceuticals | Paints | Cosmetics | Waste | Flavours | Inks | Waxes | Adhesives | Oils - Lubricating, Hydraulic, Base, Mineral, Used/Cooking Oils | Pastes

## Applications

<b>QC</b>	
Quality Control	Transport & Storage Regulations
CLP Regulations	Waste Disposal Regulations

Application:	Test Method:	Who Says So:
Transport Regulation	Small Scale, other Closed Cups	UN GHS; IATA; ADR; IMDG; CLP; DOT CFR 49-173.120
Aviation Turbine Fuel	Small Scale, Abel, Tag	ASTM D1655; Def Stan 91-91; IATA Guidance Material; ATA 103; AFQRJOS
Gas Turbine Fuel	Small Scale, Pensky-Martens	ASTM 2880
Diesel Fuel	Small Scale, Pensky-Martens	ASTM D975; D7467
Kerosines	Small Scale, Tag	ASTM D3699
Biodiesel	Small Scale	EN14213; EN14214; ASTM D6751
Environmental	Small Scale	EPA 1020 A&B; EPA CFR40 261.21
Fuel Oil	Small Scale, Pensky-Martens A&B	ASTM D396; ISO 8217
Naphthas	Small Scale, Tag	ASTM D3734; D3735
Health & Safety	Small Scale	OSHA 29 CFR 1910.106 & 1200; CPSC CFR 16-1500.43a
Water Borne Paints	Small Scale	ISO 3679; ISO 3680
Waste Products	Small Scale	91/689/EEC

## References

"For us the Stanhope Setaflash Point Tester is a valuable piece of kit and we'd struggle without it. Our main use is to detect whether a waste is hazardous or not, as this can have knock on effects in terms of cost, safety, compliance and potential processing routes. It's very useful to have it at hand as sending samples out for off site analysis simply would not be an option at times, due to potentially lengthy turn around times."

*Nick Richardson,  
Laboratory Supervisor,  
Castle Environmental*

## Discover the entire Setaflash Range

Seta Model:	Seta Part No:	Temperature Range:	Heating/Cooling Method:
Series 3 Activecool	33200-2	10 to 130°C	Peltier Cell
Series 3 Activecool (Corrosion Resistant)	33250-2	10 to 130°C	Peltier Cell
New Series 3 'Plus' (Auto Ramp) with touchscreen	33000-2	Ambient to 300°C *(0 to 300°C with coolant module)	External Cooling module (13870-0)
New Series 3 with touchscreen	30000-2	Ambient to 300°C *(0 to 300°C with coolant module)	External Cooling module (13870-0)
Series 3 Open Cup	31000-0	Ambient to 300°C *(0 to 300°C with coolant module)	External Cooling module (13870-0)
Series 8 High Temperature	82000-0	Ambient +5 to 300°C	Ceramic Pad, Forced Air (post-test cool down)
Series 8 High Temperature (Gas)	82050-0	Ambient +5 to 300°C	Ceramic Pad, Forced Air (post-test cool down)
Series 8 Activecool	82100-0	Air: 10 to 130°C Water Assisted: -20 to 130°C	Peltier Cell
Series 8 Activecool (Gas)	81150-0	Air: 10 to 130°C Water Assisted: -20 to 130°C	Peltier Cell
Series 8 Activecool (Corrosion Resistant)	82150-0	Air: 10 to 130°C Water Assisted: -20 to 130°C	Peltier Cell
Series 8 Activecool (Corrosion Resistant, Gas)	82160-0	Air: 10 to 130°C Water Assisted: -20 to 130°C	Peltier Cell

