

# OMEGA 3 – FISH OIL SUPPLEMENT

## DIGESTION OF OMEGA 3 FISH OIL USING COLDBLOCK™ DIGESTION TECHNOLOGY

*Author: Lorie-Anne Doig*

### Introduction

This application note will focus on the digestion of Omega 3 using ColdBlock™ Digestion Technology.

### Method

Samples were digested using the following method:

- Chiller temperature was set to -5°C
- An omega 3 pill was weighed and placed into a ColdBlock™ Digestion vessel.
- 4 mL H<sub>2</sub>SO<sub>4</sub> was added and sample was digested at 100% for 5 minutes
- Cool for 2 minutes, then slowly and carefully add 20mL HNO<sub>3</sub> and digest again at 100% for 20 minutes
- Sample was cooled and diluted 50mL using 2% HNO<sub>3</sub>.

### Instrument

ColdBlock™ Digestion CB12L Technology.

### General

This procedure is specific for the sample digested and may need modification for different samples to achieve the desired result.